

EZ:faast Amino Acid Analysis of Wine (GC/MS)

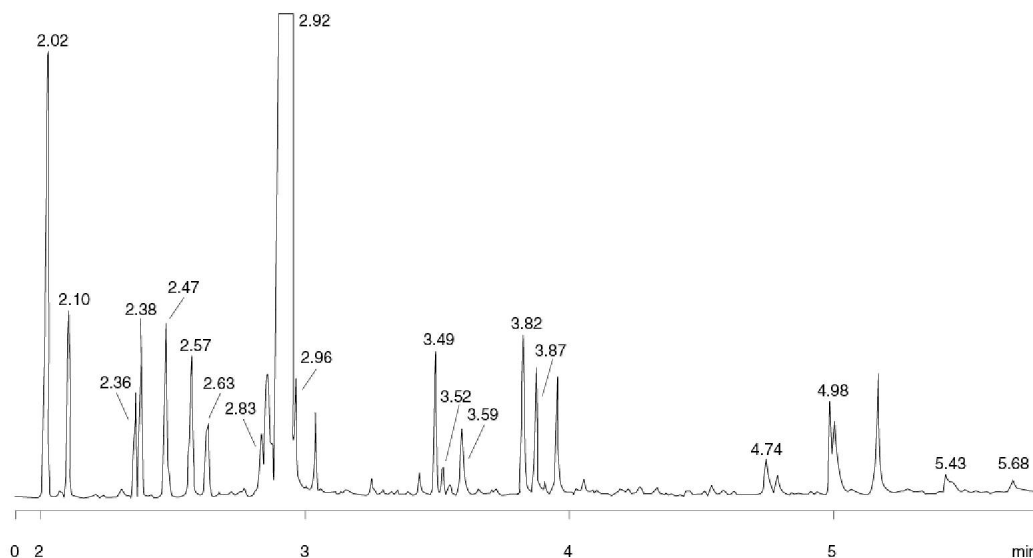
Page 1 of 1

Column: EZ:faast GC
Phase: Zebron ZB-AAA (proprietary)
Dimensions: 10 meters x 0.25 mm x 0 µm
Order No: KG0-7166
Oven Profile: 30°C/min from 110° to 320°C, hold at 320°C for 1 minute
Carrier Gas: Constant Flow Helium, 1.1 mL/min
Injection: Split 15:1 2.5 µL @ 250 °C
Detection: Mass Spectrometer (MS) (300 °C)

Products used in this Application:

EZ:faast[™]**Analyst Note:** Hot needle injection

14610

**ANALYTES:**

1 ALA	10 ASN	19 TRP
2 GLY	11 ASP	
3 VAL	12 MET	
4 BAIB	13 HYP	
5 STD	14 GLU	
6 LEU	15 PHE	
7 ILE	16 ORN	
8 THR	17 LYS	
9 PRO	18 TYR	

©Copyright 2000-2005 Phenomenex Inc.

For more information call your Phenomenex Representative at (310) 212-0555

www.phenomenex.comPhenomenex products are available worldwide. For the distributor in your country, contact Phenomenex USA, International Department by telephone, fax or e-mail: international@phenomenex.com.

USA
tel.: (310) 212-0555
fax: (310) 328-7768
email: info@phenomenex.com

Puerto Rico
(800) 541-HPLC
(310) 328-7768
info@phenomenex.com

Canada
(800) 543-3681
(310) 328-7768
info@phenomenex.com

France
01 30 09 21 10
01 30 09 21 11
franceinfo@phenomenex.com

United Kingdom
01625-501367
01625-501796
ukinfo@phenomenex.com

Germany
06021-58830-0
06021-58830-11
anfrage@phenomenex.com

New Zealand
09-4780951
09-4780952
info@phenomenex.co.nz

Australia
02-9428-6444
02-9428-6445
info@phenomenex.com.au