


Cold Pressed Orange Oil

Column:	Zebron ZB-5ms	Gold Seal:	Phenomenex single groove	(AG0-7518)
Dimensions:	20meter x 0.18mm x 0.18µm	Liner:	Single Taper/Gooseneck	(AG0-4680)
Order No.:	7FD-G010-08	Septum:	Phenomenex PhenoRed-400	(AG0-4690)
Injection:	Split 100:1 @ 180°C, 1µL			
Carrier Gas:	Constant Flow Helium 1.0 mL/min			
Oven Program:	50°C for 2 min to 100°C at 8°C/min to 225°C at 18°C/min for 4 min			
Detector:	MSD; 30-350 amu			
Sample:	1. α-Pinene 2. β-Phellandrine 3. β-Myrcene 4. Octanal 5. 3-Carene 6. Limonene 7. Linalool 8. Nonanal 9. <i>cis</i> -Limonene oxide 10. <i>trans</i> -Limonene oxide 11. Citronellal 12. α-Terpineol 13. Decanal 14. Neral 15. Carvone 16. Geranial 17. α-Cubenene 18. β-Cubenene 19. Dodecanal 20. Valencene 21. Cadinine 22. Nootkatone			

Notes: Sample was 10% in Dichloromethane

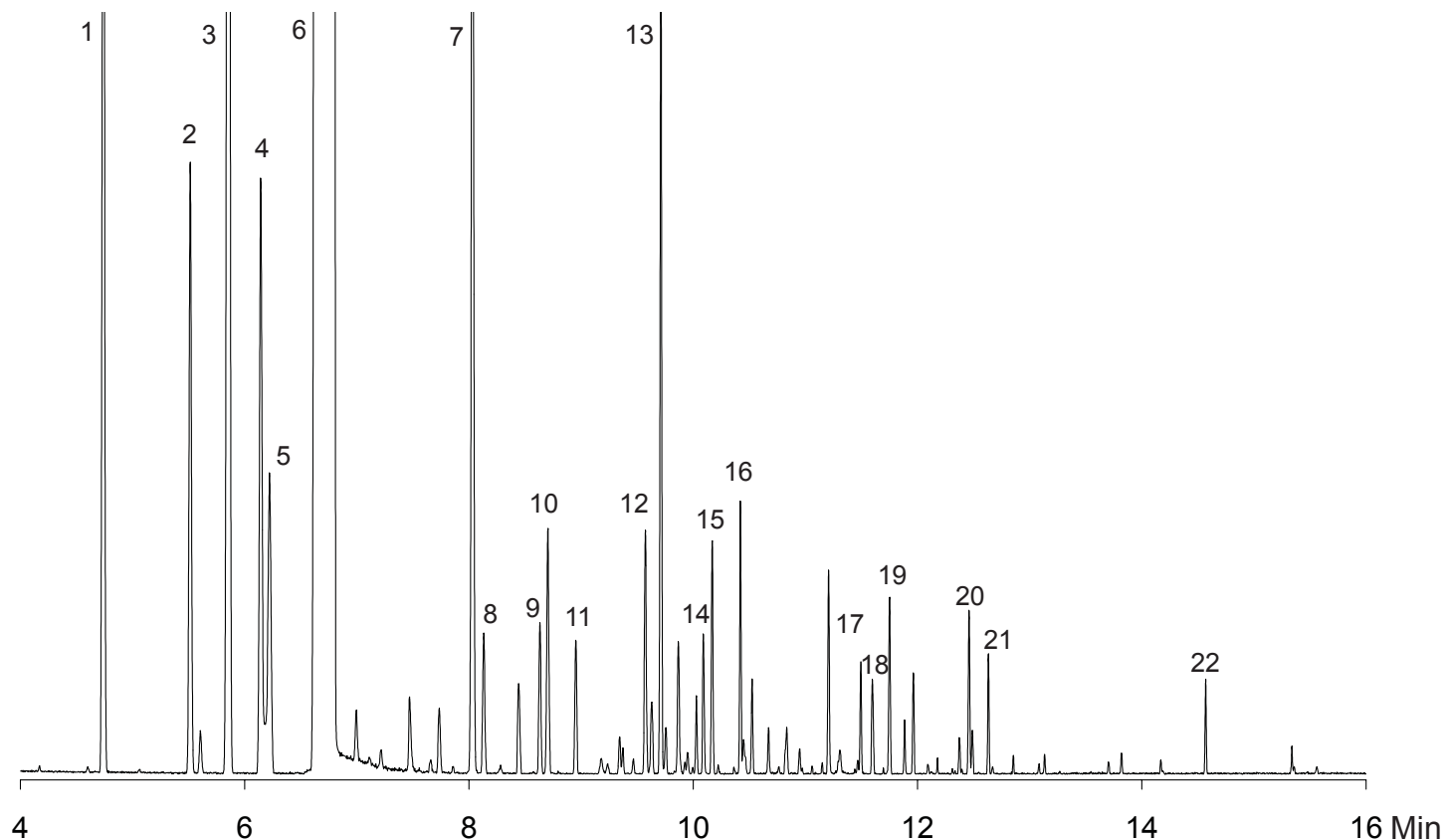

Phenomenex products are available worldwide. For the distributor in your country, contact Phenomenex USA, International Department by telephone, fax or e-mail: international@phenomenex.com.

phenomenex
...breaking with tradition™

USA	Puerto Rico	Canada	United Kingdom	Germany	New Zealand	Australia
mail: 411 Madrid Avenue Torrance, CA 90501-1430 USA	273 Sierra Morena, Suite #104 San Juan, Puerto Rico 00926	411 Madrid Avenue Torrance, CA 90501-1430 USA	Queens Avenue, Hurdfield Ind. Est., Macclesfield, Cheshire SK10 2BN, England	Zeppelinstr. 5 63741 Aschaffenburg Deutschland	P.O. Box 31-601 Milford Auckland New Zealand	P.O. Box 4084 Lane Cove, NSW 1595 Australia
tel.: (310) 212-0555 fax: (310) 328-7768 email: info@phenomenex.com	(800) 541-HPLC (310) 328-7768 info@phenomenex.com	(800) 543-3681 (310) 328-7768 info@phenomenex.com	01625-501367 01625-501796 ukinfo@phenomenex.com	06021-58830-0 06021-58830-11 anfrage@phenomenex.com	09-4780951 09-4780952 info@phenomenex.co.nz	02-9428-6444 02-9428-6445 info@phenomenex.com.au