


# SPE VACUUM MANIFOLDS

## Operating & Assembly Instructions

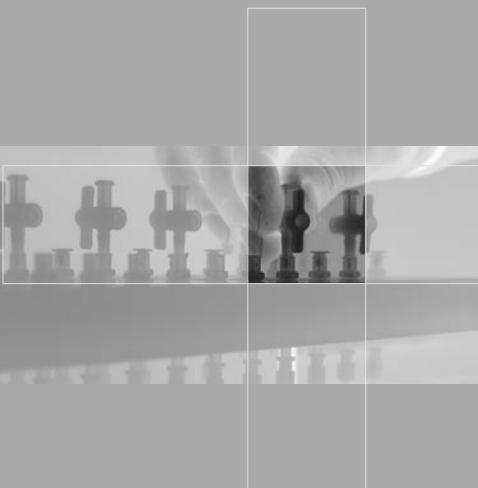

[www.phenomenex.com](http://www.phenomenex.com)  
Phenomenex products are available worldwide. For the distributor in your country,  
contact Phenomenex USA, International Department by telephone, fax or e-mail: [international@phenomenex.com](mailto:international@phenomenex.com).



...breaking with tradition<sup>™</sup>

USA		Puerto Rico		Canada		France		United Kingdom		Germany		New Zealand		Australia	
mail:	411 Madrid Avenue Torrance, CA 90501-1450 USA	275 Sierra Morena, Suite #104 San Juan, Puerto Rico 00926		411 Madrid Avenue Torrance, CA 90501-1450 USA		118130 avenue Jean Jaures 75317 Paris Cedex 19 France		Queens Avenue Hurstfield, Ess., Melbourne, Victoria 3100 20th, England		Zappelnstr. 5 65741 Haschenburg Deutschland		P.O. Box 31601 Milton Auckland New Zealand		P.O. Box 4084 Lane Cove, NSW 2066 Australia	
tel:	310-212-0555	800-541-HPLC 310-328-7768		800-543-3681 310-328-7768		01 72 72 92 10 01 72 72 92 11		01 625-501367 01 625-501796		06021-58830-0 06021-58830-11		09-4780951 09-4780952		02-9428-6444 02-9428-6445	
fax:	310-328-7768														
e-mail:	info@ phenomenex.com	info@ phenomenex.com		info@ phenomenex.com		franceinfo@ phenomenex.com		ukinfo@ phenomenex.com		airfrage@ phenomenex.com		info@ phenomenex.co.nz		info@ phenomenex.com.au	

# NEW

## SPE Tubes From Phenomenex



# strata™

**Tired of clogging and irreproducible SPE tubes? Try Strata™ solid phase extraction products manufactured by Phenomenex.**

strata™ features an extremely narrow particle size range. When particle size distribution is well controlled, your sample and solvent will flow uniformly through the sorbent bed, maximizing sample-sorbent contact. The benefits are optimum flow rates and high recoveries. Tube-to-tube flow is also more consistent, making sample processing on vacuum manifolds and automated devices more reliable and less prone to clogging or drying out.

Over 25 different tests are performed on each batch of strata including tests for: purity, recovery, and flow. QC data obtained throughout the production process is documented to provide complete traceability. You can be confident the strata products you order 6 months, 1 year or 2 years from now will perform just like the strata products you order today.

### **Product Selection & Availability**

strata tubes are constructed of ultra-pure medical grade polypropylene, Luer-tipped syringe barrels with 1, 3, and 6mL reservoir volumes. Sorbent chemistries include a complete range of high performance non-polar, polar and ionic phases, both silica and polymer based in masses ranging from 30 to 1000mg.

strata products are available through all Phenomenex offices and authorized distributors, for fast delivery anywhere in the world.

**Please call for ordering information.**

# SPE

## Reference Manual & Users Guide

New from Phenomenex, an invaluable tool for the SPE novice and expert alike. Includes:

- Protocols for reversed, normal phase and ion exchange extractions
- Method development & optimization tips
- Sorbent selection flowchart



**Hard copy version**  
**Order No.: AAO-6067**  
62 pages in length


Order No.	Manifold Set	Unit
AH0-6048	Stopcocks, for 12-position manifold	12/pk
AH0-6049	Stopcocks, for 24-position manifold	24/pk
AH0-6050	Drying Attachment, for 12-position manifold	ea
AH0-6051	Drying Attachment, for 24-position manifold	ea
AH0-6052	12-Position Vacuum Waste Container, polypropylene	10/pk
AH0-6053	Accessory, Female Luer Fittings	2/pk
AH0-6054	Accessory, Male Luer Fittings	2/pk
AH0-6055	Accessory, Support Posts for Rack	3/pk
AH0-6056	Accessory, Legs for Cover, black	4/pk
AH0-6057	Accessory, Vacuum Gauge and Valve Assembly	ea
AH0-6058	Accessory, Valve Assembly Only	ea
AH0-6059	Accessory, Vacuum Gauge Only	ea
AH0-6060	Accessory, Retaining Clips	12/pk
AH0-6061	Accessory, Vacuum Manifold Plugs	50/pk
AH0-6062	Accessory, Control Valve	25/pk
AH0-6063	Accessory, Control Valve	50/pk
AH0-6064	Accessory, Teflon Needles 100/pk	
AH0-6065	Accessory, Teflon Needles 500/pk	
AH0-7191	Accessory, Adaptor Caps for 1, 3 and 6mL SPE tubes	15/pk
AH0-7378	Accessory, Adaptor Caps for 12 and 20mL SPE tubes	5/pk

- NOTES:**
1. The 10-position Tall Boy Vacuum Manifold Collection Rack includes 4 plates: one base plate, one dimple plate, one small plate and one large plate and three riser legs, along with 12 manifold clips to support the plates. The assembly also includes 10 polypropylene needles, 10 stopcocks and 4 black legs to support the lid when taken off the glass block.
  2. The 12-position Collection Rack consists of 3 support posts, bottom plate, 13mm plate, 16mm plate, autosampler plate, volumetric plate, and 12 retaining clips.
  3. The 24-position Collection Rack consists of 3 support posts, bottom plate, dimple plate, 13mm plate, 16mm plate, and

# SPE VACUUM MANIFOLDS

## Operating & Assembly Instructions

**CAUTION:** Do not operate this manifold without first reading and fully understanding the operating instructions.

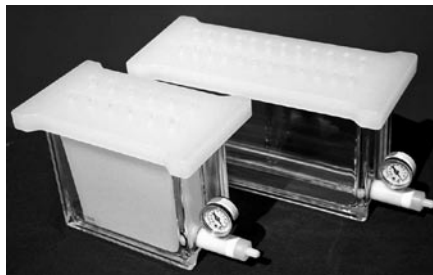

## INTRODUCTION

Phenomenex SPE vacuum manifolds are available in 10-, 12- and 24-position configurations. These manifolds permit convenient batch processing of SPE or sample filtration and help ensure consistent extraction and filtration results. The manifolds consist of a clear glass chamber and lid to which a vacuum is applied thereby drawing the sample through the SPE cartridge or filtration device. Adjustable racks placed in the glass vacuum chamber accommodate a variety of sample collection vessels and test tube configurations.

## Table of Contents

Manifold Assembly .....	2
Vacuum Source and Waste Trap .....	4
Manifold Operation .....	5
Optional Accessories .....	6



figure 1

1. Lid Assembly:

A. Attach the four black legs to the underside of the manifold lid as shown in figure 1.



figure 2

B. Attach the needles to the male luer connection fittings on the underside of the lid as shown in figure 2.



figure 3

C. Firmly insert the stopcocks into the female luer fittings on top of the lid as shown in figure 3.

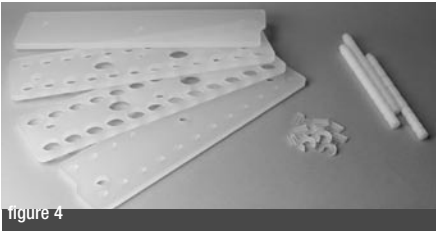


figure 4

2. Rack Assembly:

The rack assembly consists of a base, three attachment posts, a dimpled tube support platform and an upper rack with either 16mm or 12mm diameter holes. Select the upper rack with

Order No.	Replacement Parts	Unit
AH0-6027	Cover, Gasket and 12 Stopcocks for 12-position manifold	ea
AH0-6028	Cover, Gasket and 24 Stopcocks for 24-position manifold	ea
AH0-6029	Gaskets for 12-position manifold	2/pk
AH0-6030	Gaskets for 24-position manifold	2/pk
AH0-6031	Vacuum Gauge, Valve and Glass Chamber for 12-position manifold	ea
AH0-6032	Vacuum Gauge, Valve and Glass Chamber for 24-position manifold	ea
AH0-6033	Manifold Needles, polypropylene	12/pk
AH0-6034	Manifold Needles, polypropylene	24/pk
AH0-6035	Manifold Needles, stainless steel	12/pk
AH0-6036	Manifold Needles, stainless steel	24/pk
AH0-6037	Collection Rack, including shelves, legs, clips and posts, for 12-position manifold	ea
AH0-6038	Collection Rack, including shelves, legs, clips and posts, for 24-position manifold	ea
AH0-6040	Plate for 13mm Test Tube, for 12-position manifold	ea
AH0-6041	Plate for 13mm Test Tube, for 24-position manifold	ea
AH0-6042	Plate for Volumetric Flask, for 12-position manifold	ea
AH0-6043	Plate for 16mm Test Tube, for 12-position manifold	ea
AH0-6044	Plate for 16mm Test Tube, for 24-position manifold	ea
AH0-6045	Plate for Autosampler Vial, for 12-position manifold	ea
AH0-6046	Plate, Dimple, for 12-position manifold	ea
AH0-6047	Plate, Base, for 12-position manifold	ea

**1. Teflon® Needles and Teflon® Control Valves**  
 incorporating flow control are available. These are recommended for trace analysis requiring a plasticizer-free, inert fluid path.

**2. Drying Attachment:**  
 Optional drying attachments for the 12- and 24-position manifolds offer a convenient and efficient way to evaporate and further concentrate the SPE extracts. In addition, the drying attachment may be used to apply a positive pressure to the tops of the cartridges which is useful for drying the cartridges during the extraction procedure.

**3. Waste Container:**  
 (12-position manifold only) An optional solvent-resistant polypropylene liner is available for the 12-position manifold which can collect solvent wastes generated during the SPE extraction. Using the waste liner can greatly simplify manifold clean-up and maintenance.

**ORDERING INFORMATION**

Order No.	Manifold Set	Unit
AH0-7502	10-Position Tall Boy Vacuum Manifold, complete assembly <sup>1</sup>	ea
AH0-6023	12-Position Vacuum Manifold Set, complete assembly <sup>2</sup>	ea
AH0-6024	24-Position Vacuum Manifold Set, complete assembly <sup>3</sup>	ea
<b>Replacement Parts</b>		
AH0-7503	10-Position Glass Chamber	ea
AH0-6025	12-Position Glass Chamber	ea
AH0-6026	24-Position Glass Chamber	ea
AH0-7504	Cover, Gasket and 10 Stopcocks for 10-position Tall Boy manifold	ea

holes to match collection tube diameter. Vertical adjustment of the dimpled support platform upon which the bottoms of the test tubes rest allows the user to configure the manifold for a



figure 5

A. Lay the base onto a flat surface oriented so that the indentation is on the bottom right. Attach the three attachment posts to the base (figure 5). Correct orientation of the indentation is necessary for the rack to clear the vacuum gauge assembly when placed inside the tank.



figure 6

B. Slide the dimpled support plate onto the posts by aligning the three holes in the plate with attachment posts. The indentation on the side of the support plate must match the indentation of the base (figure 6).

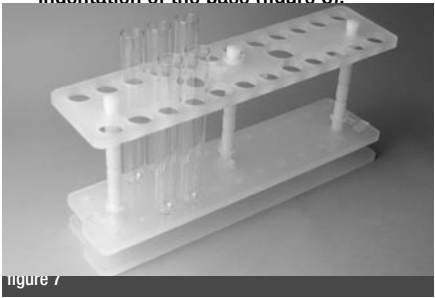


figure 7

C. Place the upper rack with the proper diameter holes onto the attachment posts as shown in figure 7.

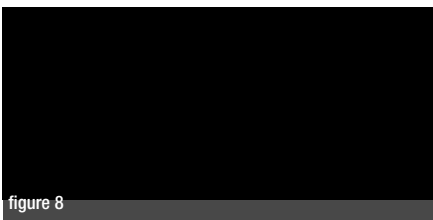


figure 8



figure 9

D. Adjust the height of the dimpled support platform shelf so that the delivery needles in the manifold lid are a minimum of 5mm inside the collection vessels (figure 8). “C”-shaped support clips attach to the posts directly under the shelf allowing height adjustment (figure 9).

E. To collect eluents, place rack loaded with test tubes into glass basin and carefully replace the lid.

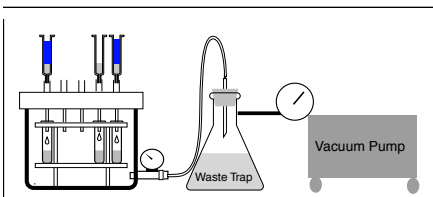


figure 10

### VACUUM SOURCE AND WASTE TRAP

Installation of a liquid waste trap between the manifold and the vacuum source is highly recommended for the protection of the vacuum source and for collection and proper disposal of extraction waste solvents. (figure 10). Use  $\frac{5}{16}$ " (8mm) ID thick wall vacuum tubing for connections between the pump, waste trap, and manifold. A high capacity, corrosion-resistant vacuum pump capable of generating a minimum vacuum of 10 inches of mercury is recommended.

### MANIFOLD OPERATION

**Caution:** Vacuum should not exceed 20 inches of Mercury (68 kPa). Exceeding this vacuum will void the manifold warranty and may cause dangerous glass basin failure.



figure 11

A. Turn on vacuum to the system and using the knurled bleed valve ring located adjacent to the manifold vacuum gauge (figure 11) adjust the maximum vacuum level. The bleed valve ring also permits rapid vacuum release of the entire system. Most analysts find that 3 to 5 inches of vacuum is satisfactory for proper flow through the SPE columns.



figure 12

by stopcocks or optional Teflon control valves. Stopcocks are fully open when the valve is in the vertical position and fully closed when in the horizontal position.

C. Individual stopcocks should be in the closed position prior to removal of the SPE column when under vacuum to prevent splashing of solvents.

D. Prior to removing the lid, it is important to completely bleed the vacuum from the manifold by opening the bleed valve ring (figure 11) or fully opening any unused stopcocks on the lid. Failure to do so may result in eluent splashing and sample cross contamination.

### OPTIONAL ACCESSORIES