




Aromatic Amines

Column: ZB-5
Dimensions: 30m x 0.25mm x 0.25µm
Flow: Constant flow of 1.0 mL/min
Injector: 220°C; 1 µL at 15:1 split
Oven: 60°C for 1min to 130°C at 12°C/min to 220°C at 30°C/min for 4min.
Detector: Full Scan MS
Analytes: 1.58 mg/mL each

Aromatic Amines on ZB-5, 30m x 0.25mm x 0.25µm

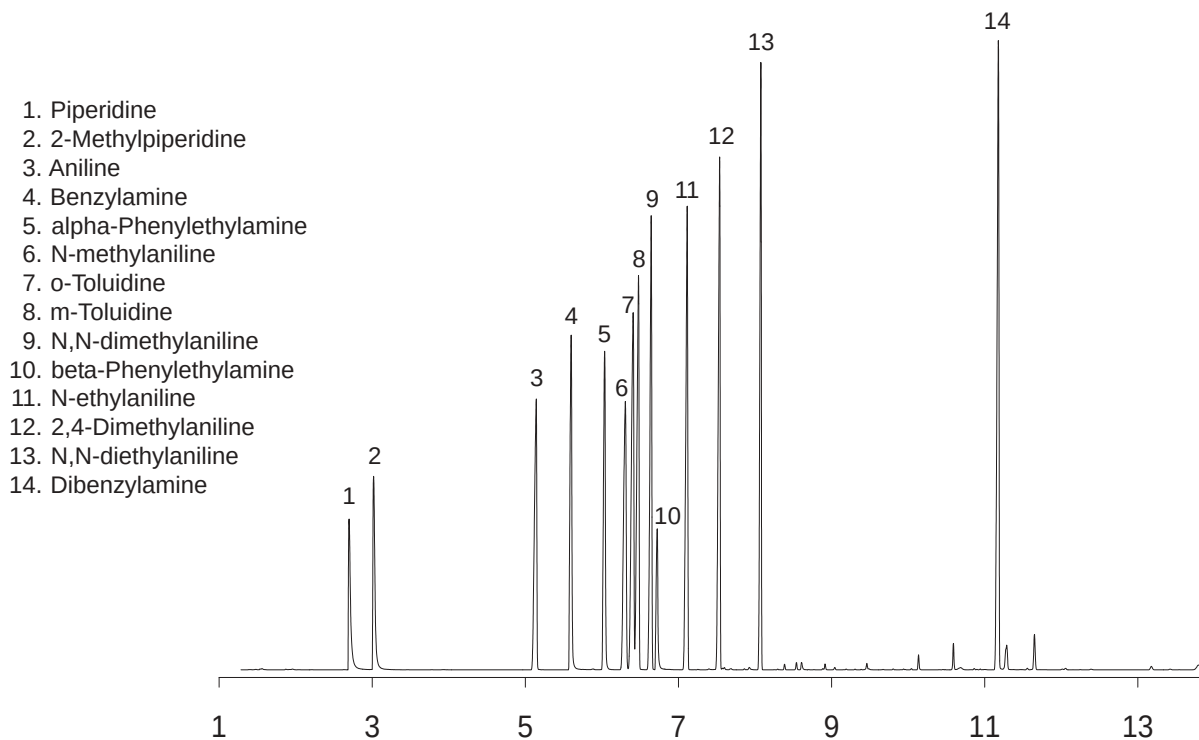

www.phenomenex.com

Phenomenex products are available worldwide. For the distributor in your country, contact Phenomenex USA, International Department by telephone, fax or e-mail: international@phenomenex.com.

phenomenex
 ...breaking with tradition


USA	Puerto Rico	Canada	United Kingdom	Germany	New Zealand	Australia
tel.: (310) 212-0555	(800) 541-HPLC	(800) 543-3681	01625-501367	06021-58830-0	09-4780951	1800-553-929
fax: (310) 328-7768	(310) 328-7768	(310) 328-7768	01625-501796	06021-58830-11	09-4780952	1800-553-923
email: info@phenomenex.com	info@phenomenex.com	info@phenomenex.com	ukinfo@phenomenex.com	anfrage@phenomenex.com	info@phenomenex.co.nz	info@phenomenex.com.au

Aromatic Amines

Column: ZB-1701
Dimensions: 30m x 0.25mm x 0.25µm
Flow: Constant flow of 1.0 mL/min
Injector: 220°C; 1 µL at 15:1 split
Oven: 60°C for 1min to 110°C at 30°C/min to 135°C at 9°C/min to 260°C at 30°C/min for 2min.
Detector: Full Scan MS
Analytes: 1.58 mg/mL each

Aromatic Amines on ZB-1701, 30m x 0.25mm x 0.25µm

1. Piperidine
2. 2-Methylpiperidine
3. Aniline
4. Benzylamine
5. alpha-Phenylethylamine
6. N-methylaniline
7. o-Toluidine
8. m-Toluidine
9. N,N-dimethylaniline
10. beta-Phenylethylamine
11. N-ethylaniline
12. 2,4-Dimethylaniline
13. N,N-diethylaniline
14. Dibenzylamine

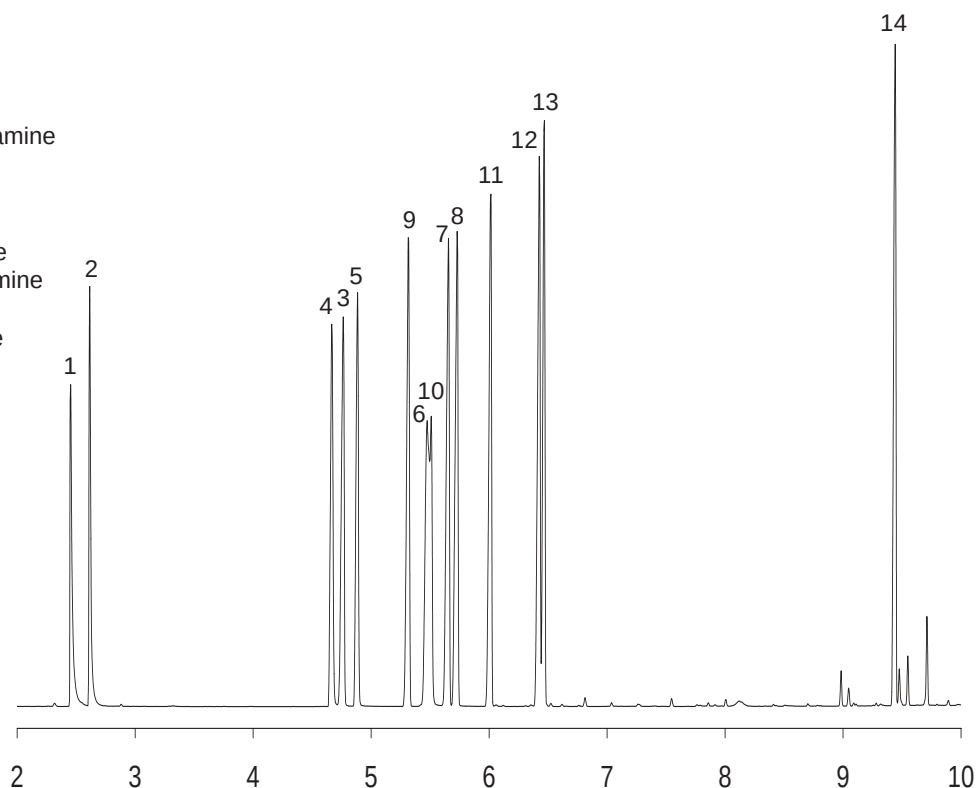

Ordering Information

Order Number	Description
7HG-G002-11	ZB-5 - 30m x 0.25mm x 0.25µm
7HG-G006-11	ZB-1701 - 30m x 0.25mm x 0.25µm