

Reported by User: System


Project Name: Vitamin

# SAMPLE INFORMATION

Sample Name: V\_B group  
Sample Type: Standard  
Vial: 1  
Injection #: 1  
Injection Volume: 20.00 ul  
Run Time: 35.0 Minutes  
Sample Set Name lee

Acquired By: System  
Date Acquired: 2006-09-21 오후 6:39:36  
Acq. Method Set: V\_B group  
Date Processed: 2006-09-22 오전 9:25:51  
Processing Method V\_B Group  
Channel Name: 2487Channel 1  
Proc. Chnl. Descr.:

## Auto-Scaled Chromatogram

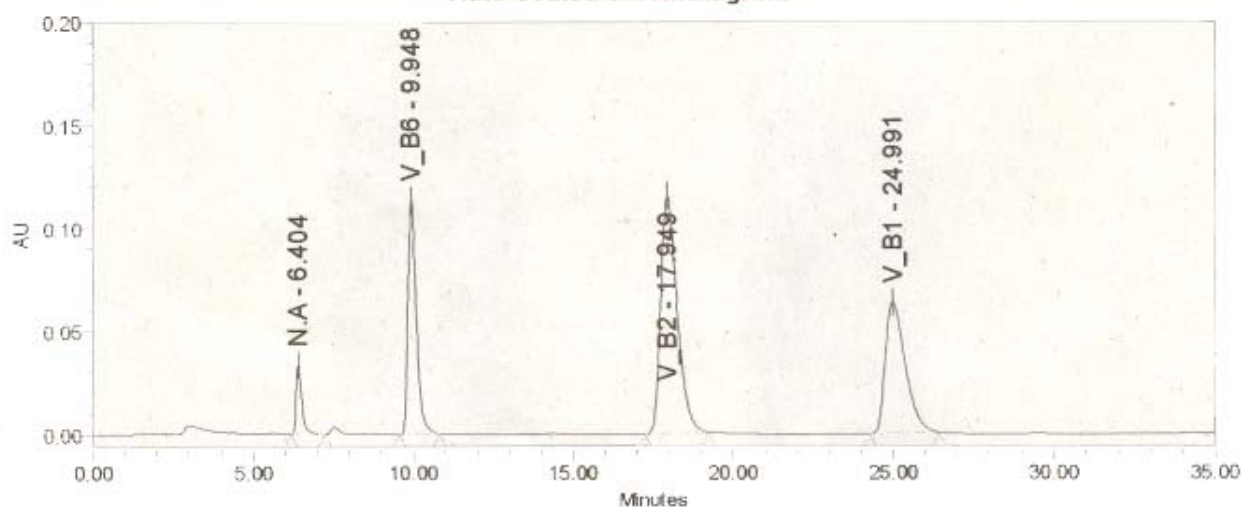

## Peak Results

|   | Name | RT     | Area    | Height | Amount | Units |
|---|------|--------|---------|--------|--------|-------|
| 1 | N.A  | 6.404  | 416746  | 33607  |        |       |
| 2 | V_B6 | 9.948  | 2072602 | 113405 |        |       |
| 3 | V_B2 | 17.949 | 4091935 | 115033 |        |       |
| 4 | V_B1 | 24.991 | 2745211 | 63496  |        |       |

M.P: 25% MeOH + B6 1g + 2% HoAc

UV: 280nm

flow: 0.6ml/min