


SPE Drying Attachment

Used with the following Part Number(s):

Catalog_No	Description_1	Description_2	UOM
AH0-6023	SPE 12-Position Vacuum	Manifold Set - Complete	Ea
AH0-6024	SPE 24-Position Vacuum	Manifold Set - Complete	Ea

Instructions:

AH0-6050 Drying Attachment

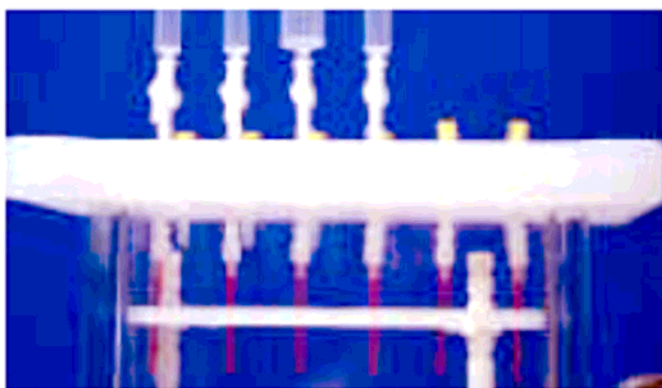

The drying attachment is used in conjunction with 12, 16 and 24 Port Manifolds to concentrate samples that have been eluted from Solid Phase Extraction columns. They can also be used to dry packing beds in SPE columns prior to elution of the sample.

Instructions To Dry Samples:

After the samples are collected in the receiving tubes inside the vacuum manifold, remove the lid and replace it with the drying attachment. Connect the drying attachment to an inert gas source by securing the tubing from the vacuum to the brass hose barb on the drying attachment. Gas then flows into each receiving tube to dry or concentrate the samples.

Instructions To Dry Packing Beds:

Leave the SPE tubes on the vacuum manifold lid. Attach an adaptor to each tube to prevent outside air from entering the tube. Attach the drying attachment lid to the adaptors on the tube with the male luers of the drying attachment. The drying attachment will make a tight connection with the column adaptor by way of the male luers. Attach an inert gas source to the drying attachment by attaching the tubing from the vacuum to the brass hose barb on the drying attaching. By applying vacuum to the manifold, the inert gas will be pulled through the tubes and dry the packing beds prior to elution of your samples.



www.phenomenex.com

Phenomenex products are available worldwide. For the distributor in your country, contact Phenomenex USA, International Department by telephone, fax or e-mail: international@phenomenex.com.

USA	Puerto Rico	Canada	United Kingdom	Germany	New Zealand	Australia
mail: 411 Madrid Ave. Torrance, CA 90501-1430 USA tel.: (310) 212-0555 fax: (310) 328-7768 email: info@phenomenex.com	271 Sierra Morena, Suite #104 San Juan, Puerto Rico 00926 (800) 541-HPLC (310) 328-7768 info@phenomenex.com	411 Madrid Ave. Torrance, CA 90501-1430 USA (800) 543-3681 (310) 328-7768 info@phenomenex.com	Queens Avenue, Hurdfield Ind. Est., Macclesfield, Cheshire SK10 2BN, England 01625-501367 01625-501796 ukinfo@phenomenex.com	Zeppelinstr. 5 63741 Aschaffenburg Deutschland 06021-58830-0 06021-58830-11 anfrage@phenomenex.com	P.O. Box 31-601 Milford Auckland New Zealand 09-4780951 09-4780952 info@phenomenex.co.nz	P.O. Box 264 Pennant Hills NSW 1715 Australia 1800-553-929 1800-553-923 info@phenomenex.com.au