

How to Convert from Packed to Capillary GC Columns!

Sky Countryman, GC Product Manager
Phenomenex Inc., Torrance, CA, USA

If you're still wondering why you should convert to a high efficiency capillary GC system, consider these three reasons:


1 EASY . . .
Takes less than 30min!

2 AFFORDABLE . . .
Costs less than \$100


3 BETTER RESULTS . . .
Faster, better resolution, improved results.

See how simple it is!


Packed Inlet

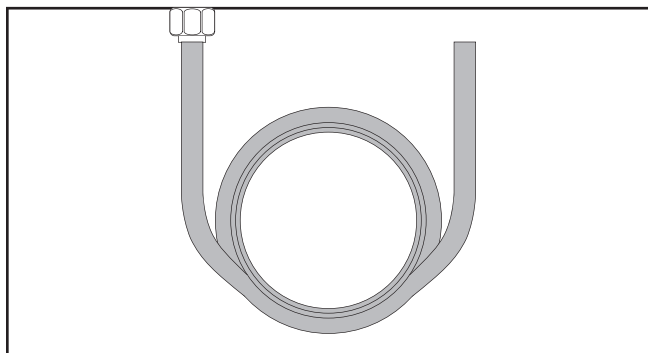


Capillary Inlet*

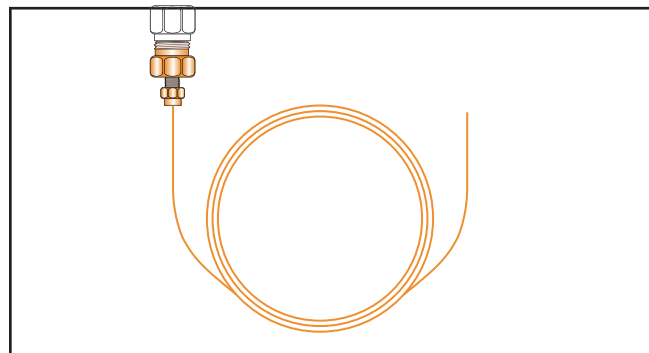

Packed Detector

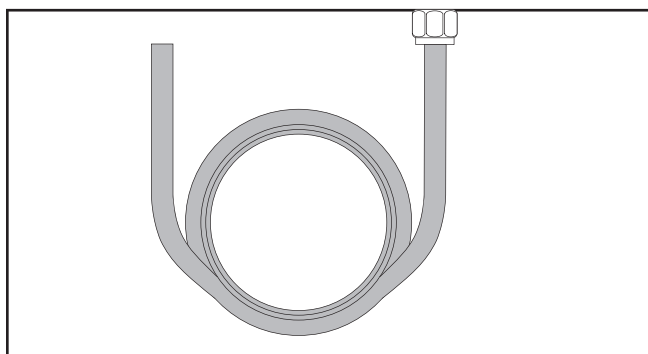

Capillary Detector*

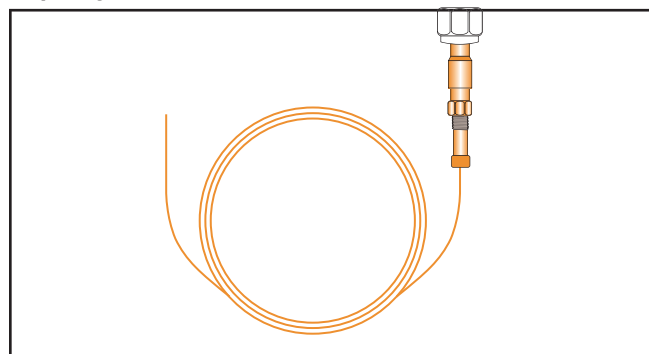

* Parts necessary for capillary conversion are those shown in orange. Conversions kits are available for all major GC brands.



www.phenomenex.com

Phenomenex products are available worldwide. For the distributor in your country, contact Phenomenex USA, International Department by telephone, fax or e-mail: international@phenomenex.com.

phenomenex
...breaking with traditionSM

USA
tel.: (310) 212-0555
fax: (310) 328-7768
email: info@phenomenex.com

Puerto Rico
(800) 541-HPLC
(310) 328-7768
info@phenomenex.com

Canada
(800) 543-3681
(310) 328-7768
info@phenomenex.com

France
01 30 09 21 10
01 30 09 21 11
franceinfo@phenomenex.com

United Kingdom
01625-501367
01625-501796
ukinfo@phenomenex.com

Germany
06021-58830-0
06021-58830-11
anfrage@phenomenex.com

New Zealand
09-4780951
09-4780952
info@phenomenex.co.nz

Australia
02-9428-6444
02-9428-6445
info@phenomenex.com.au

Benefits of Capillary Vs. Packed Columns

1. Better Resolution:


- Packed GC ~ 1,800-3,600 plates/column
- Capillary GC ~ 58,500 - 234,000 plates/column

2. Faster Analysis

3. Lower Bleed


4. Better Reproducibility

2meter x 2mm ID Glass Packed Column

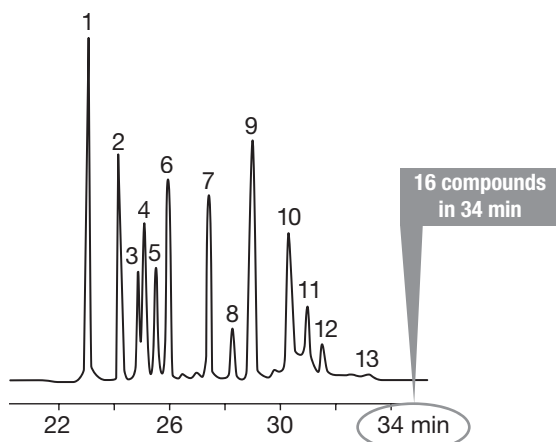

1. α -BHC
2. γ -BHC
3. β -BHC
4. Heptachlor
5. δ -BHC
6. Aldrin
7. Heptachlorepoxide
8. Endosulfan I
9. Dieldrin/4,4'-DDE
10. Endrin
11. Endosulfan II/4,4'-DDD
12. 4,4'-DDT
13. Endosulfan sulfate/Endrin aldehyde

30 meter x 0.25mm ID x 0.25 μ m Capillary Column APP ID 15785

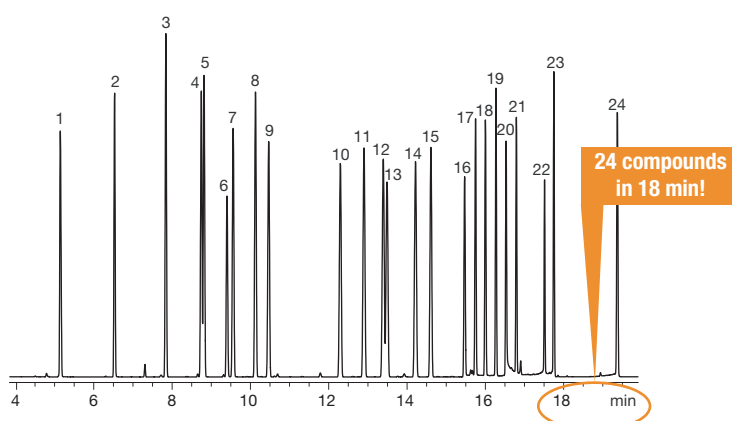

1. 1-Bromo-2-Nitrobenzene (IS)
2. Tetrachlorometaxylene (surr)
3. α -BHC
4. Pentachloronitrobenzene (surr)
5. γ -BHC
6. β -BHC
7. Heptachlor
8. δ -BHC
9. Aldrin
10. Heptachlor Epoxide
11. γ -Chlordane
12. α -Chlordane
13. Endosulfan I
14. DDE
15. Dieldrin
16. Endrin
17. DDD
18. Endosulfan II
19. DDT
20. Endrin Aldehyde
21. Endosulfan Sulfate
22. Methoxychlor
23. Endrin Ketone
24. Decachlorobiphenyl (IS)

For more information about the benefits of using capillary GC columns or to find out what you need to convert your system to use capillary columns, please contact your local Phenomenex representative.