


APPLICATION NOTE - HPLC
Estradiol - USP Method



15001

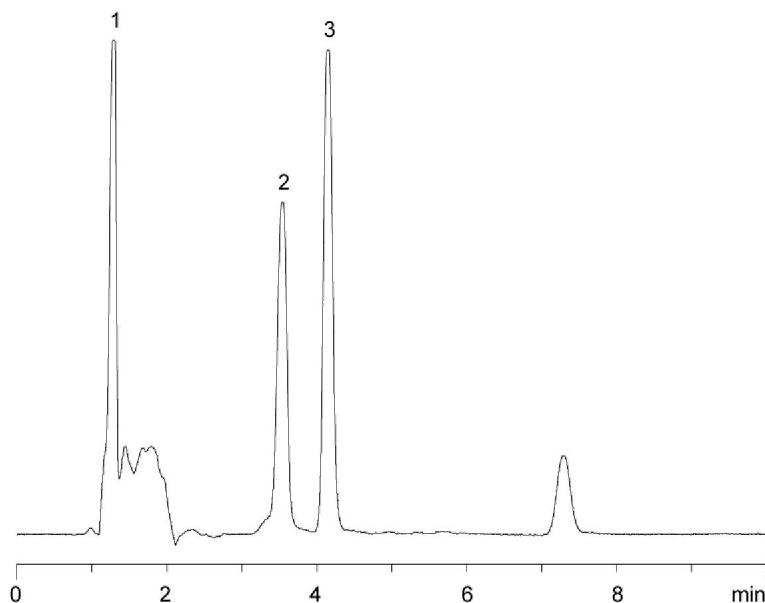

Column: Gemini 5 μ m C18 / ODS C18

Dimensions: 150 x 4.6 mm ID

Order No: 00F-4435-E0

Elution Type: Isocratic

Eluent A: Water

Pct. A: 45

Eluent B: Acetonitrile

Pct. B: 55

Flow Rate: 1 mL/min

Temperature: 25 °C

Detection: UV-Vis Abs. - Variable Wave. @ 205 nm

SAMPLE ANALYTES:

- 1 Estrone
- 2 Ethyl paraben
- 3 Estradiol

SPE:

For more information call your Phenomenex Representative at

©Copyright 2000-2005 Phenomenex Inc.

Ref. No.: 15001/BRO

Page 1 of 1

