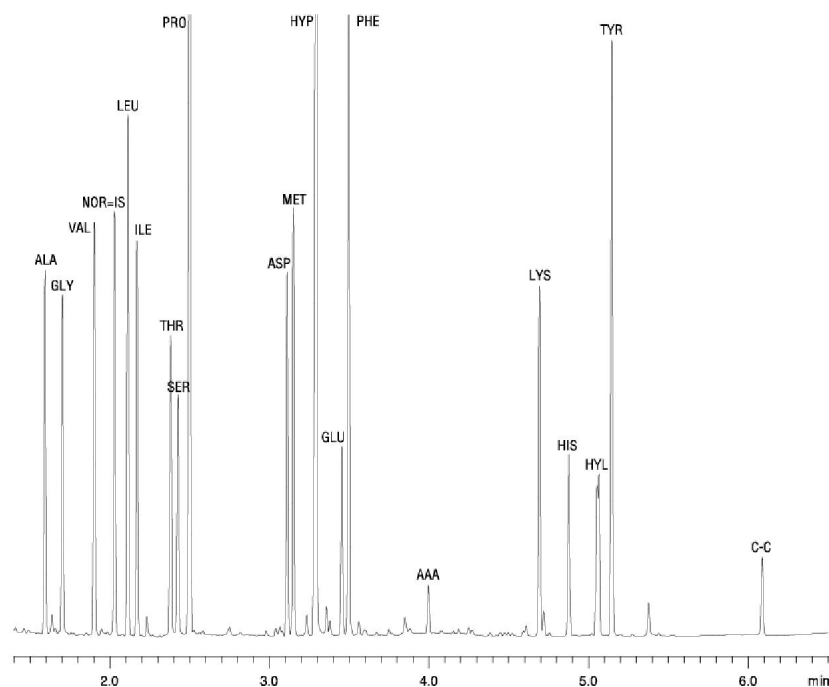


Protein Amino Acid Standards

Page 1 of 1

Column: EZ:faast GC
Phase: Zebron ZB-AAA (proprietary)
Dimensions: 10 meters x 0.25 mm x 0 µm
Order No: KG0-7167
Oven Profile: 30°C/min from 110° to 320°C, hold at 320°C for 1 minute
Carrier Gas: Constant Flow Helium, 1.47 mL/min
Injection: Split 15:1 2 µL @ 250 °C
Detection: Flame Ionization (FID) (320 °C)

Products used in this Application:

EZ:faast™


ANALYTES:

6 ALA	15 ASP	24 C-C
7 GLY	16 MET	
8 VAL	17 HYP	
9 INTERNAL STD	18 GLU	
10 LEU	19 PHE	
11 ILE	20 LYS	
12 THR	21 HIS	
13 SER	22 HYL	
14 PRO	23 TYR	

©Copyright 2000-2005 Phenomenex Inc.

For more information call your Phenomenex Representative at (310) 212-0555



www.phenomenex.com
 Phenomenex products are available worldwide. For the distributor in your country,
 contact Phenomenex USA, International Department by telephone, fax or e-mail: international@phenomenex.com.


 ...breaking with tradition™

USA tel.: (310) 212-0555 fax: (310) 328-7768 email: info@phenomenex.com	Puerto Rico (800) 541-HPLC (310) 328-7768 info@phenomenex.com	Canada (800) 543-3681 (310) 328-7768 info@phenomenex.com	France 01 30 09 21 10 01 30 09 21 11 franceinfo@phenomenex.com	United Kingdom 01625-501367 01625-501796 ukinfo@phenomenex.com	Germany 06021-58830-0 06021-58830-11 anfrage@phenomenex.com	New Zealand 09-4780951 09-4780952 info@phenomenex.co.nz	Australia 02-9428-6444 02-9428-6445 info@phenomenex.com.au
---	---	--	--	--	---	---	--